

# SARS-CoV-2 (Spike Protein) Peptide Pool

SARS-CoV-2 (spike protein) peptide pool for immune cell activation; contains two subpools that span the full-length spike protein

Catalog #100-0676

2 Vials

~25 µg (15 nmol)/peptide

## Product Description

SARS-CoV-2 (Spike Protein) Peptide Pool is provided as two lyophilized mixtures (subpools) from the SARS-CoV-2 spike protein. Each subpool contains 158 peptides, for a total of 316 peptides. The virus attaches to the cell membrane of the host through the interaction between spike protein and angiotensin-converting enzyme 2 (ACE2) receptor, and the spike protein plays a critical role in viral entry (Hoffmann et al.; Walls et al.). The subpools consist of 15-mer peptides with 11-amino-acid overlaps that cover amino acids 1 - 1273 on the spike protein. One unit of this product (i.e. ~25 µg/peptide) is sufficient for stimulating  $2.5 \times 10^8$  cells.

## Product Information

**Amino Acid Sequence:** MFVFLVLLPLVSSQCVNLTTRTQLPPAYTNSFTRGVYYPDKVFRSSVLHSTQDLFLPFFSNVTWFHAIHVSGT  
NGTKRFDNPLPFNDGVYFASTSEKSNIRGWIFGTTLDSTKQSLIVNNATNVVIVKVEFCQFCNDPFLGVYVYH  
KNNKSWMESEFRVYSSANNCTFEYVSQPFLMDLEGKQGNFKNLREFVFNKIDGYFKIYKHTPINLVRDLP  
QGFSALEPLVDLPIGINITRFQTLALHRSYLTGPDSSSGWTAGAAAYVGYLQPRFTLLKYNNENGTITDAVDC  
ALDPLSEKCTLKSFTVEKGIYQTSNFRVQPTESIVRFPNITNLCPFGEVFNATRFASVYAWNRKRISNCVADY  
SVLYNSASFSTFKCYGVSPTKLNLDLCTNRYADSFVIRGDEVRQIAPGQTGKIADYNYKLPDDFTGCVIAWNS  
NNLDSKVGGNYNLYRFRKSNLKPFERDISTEIQAGSTPCNGVEGFNCYFPLQSYGFQPTNGVGYQPYRV  
VVLSFELLHAPATVCGPKKSTNLVKNKCVNFNFNGLTGTGVLTESNKKFLPFQFGRDIADTTDAVRDPQTL  
EILDITPCSFGGVSVITPGTNTSNQVAVLYQDVNCTEVPVAIHADQLTPTWRVYSTGNSVVFQTRAGCLIGAEHVNNYSY  
ECDIPGAGICASYQTQTSNPRRARSVASQSIIAYTMSLGAENSVAYSNNNSIAIPTNFTISVTTEILPVSMTKTSV  
DCTMYICGDSTECSNLLLQYGSFCTQLNRALTGIAVEQDKNTQEVFAQVKQIYKTPPIKDFGGFNFSQILPDP  
SKPSKRSFIEDLLFNKVTLADAGFIKQYGDCLGDIARDLCAQKFNGTLVPLLLTDEMIAQYTSALLAGTITS  
GWTFGAGAALQIPFAMQMAYRFNGIGVTQNVLYENQKLIANQFNSAIGKIQDLSSTASALGKLQDVVNQN  
AQALNTLVKQLSSNFGAISSVNDILSRLDKVEAEVQIDRLITGRQLSLQTYVTQQQLIRAAEIRASANLAATKMS  
ECVLGQSKRVDFCGKGYHLMSPQSAHPGVVFLHVTYVPAQEKNFHTTAPAICHDKAHFPREGVFSVNGT  
HWFVTQRNFYEPQIITDNTFVSGNCDVVIGVNNNTVYDPLQPELDSFKEELDKYFKNHTSPDVLGDISGI  
NASVVNIQKEIDRLNEVAKNLNESLIDLQELGKYEQYIKWPWYIWLGFIAGLIAIVMVTIMLCCMTSCCCLK  
GCCSCGSCCKFDEDDSEPVKGVKLYHT

**Product Formulation:** Lyophilized as trifluoroacetate salts

**Source:** SARS-CoV-2 (severe acute respiratory syndrome coronavirus 2)

**Number of Peptides:** 158 per vial (316 total)

## SARS-CoV-2 (Spike Protein) Peptide Pool

Protein ID:	PODTC2
Protein Name:	Spike glycoprotein, S1 protein
Gene Name:	S
Purity:	Average 70%

## Preparation and Storage

**Stability and Storage:** Store at -20°C. Stable as supplied until expiry date (EXP) on label.  
NOTE: SARS-CoV-2 (Spike Protein) Peptide Pool 1 (Catalog #300-0280) and SARS-CoV-2 (Spike Protein) Peptide Pool 2 (Catalog #300-0281) are sold as part of the SARS-CoV-2 (Spike Protein) Peptide Pool (Catalog #100-0676) and are not available for individual sale.

**Preparation:** Warm to room temperature (15 - 25°C) before reconstitution. Add pure dimethyl sulfoxide (DMSO; ~40 µL) and dilute with water to the desired concentration. **Combination of the two subpools after reconstitution is not recommended.** Final concentration of DMSO must be below 1% (v/v) to avoid toxicity in the biological system. If not used immediately, aliquot and store at -20°C. Protect from light. Avoid repeated freeze-thaw cycles.

## Related Products

For a complete list of cytokines or peptide pools, as well as related products available from STEMCELL Technologies, visit [www.stemcell.com/cytokines](http://www.stemcell.com/cytokines) or contact us at [techsupport@stemcell.com](mailto:techsupport@stemcell.com).

## References

Hoffmann M et al. (2020) SARS-CoV-2 Cell entry depends on ACE2 and TMPRSS2 and is blocked by a clinically proven protease inhibitor. Cell 181(2): 271–80.e8.

Walls AC et al. (2020) Structure, function, and antigenicity of the SARS-CoV-2 spike glycoprotein. Cell 181(2): 281–92.e6.

PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED.

Copyright © 2024 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, and Scientists Helping Scientists are trademarks of STEMCELL Technologies Canada Inc. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.