

# mTeSR™1

## Defined, Feeder-Independent hESC & hiPSC Maintenance Medium

### mTeSR™1 is the Most Widely Published Feeder-Independent hESC & hiPSC Maintenance Medium

mTeSR™1 is a standardized medium for the feeder-independent maintenance of hESCs and hiPSCs.<sup>1</sup> It is a complete, serum-free, defined formulation based on Ludwig et al.<sup>2</sup> and developed under license from the WiCell™ Research Institute. To date, mTeSR™1 has been used successfully to maintain over 50 independently derived hESC and hiPSC lines.

mTeSR™1 is designed to be used with BD Matrigel™ hESC-qualified Matrix (BD Catalog #354277) as a substrate. STEMCELL Technologies pre-qualifies each batch of BD Matrigel™ to ensure consistency, reproducibility and reliability in performance.

Developed at the:



| PRODUCT | QUANTITY    | CATALOG # |
|---------|-------------|-----------|
| mTeSR™1 | 500 mL      | 05850     |
|         | 10 x 500 mL | 05870     |
|         | 25 x 500 mL | 05875     |
|         | 1L          | 05857     |

NEW!

### VIDEO

See how leading researchers have  
achieved worldwide recognition with  
mTeSR™1

[www.stemcell.com](http://www.stemcell.com)

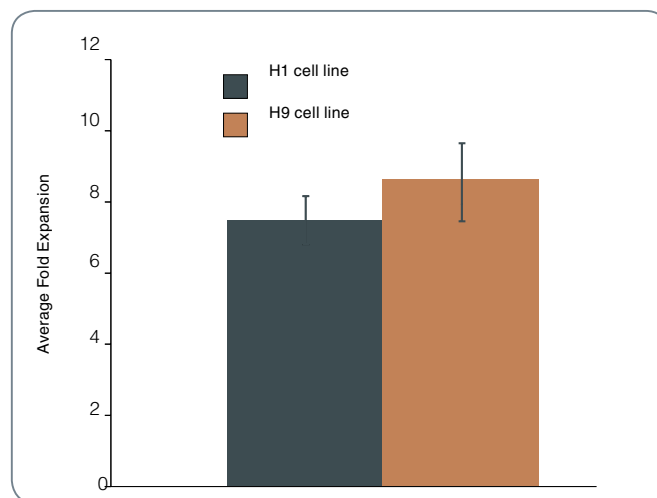
SCAN ME ▶



### Advantages of mTeSR™1

- Complete formulation that is defined and serum-free
- No feeder or conditioned media preparation
- Standardized culture method minimizes experimental variability

FIGURE 1. Consistent Expansion of hESCs & hiPSCs in mTeSR™1



H1 and H9 hESCs were expanded in mTeSR™1 for 19 and 18 passages respectively. Cultures show 7- to 10-fold expansion consistently across passages.

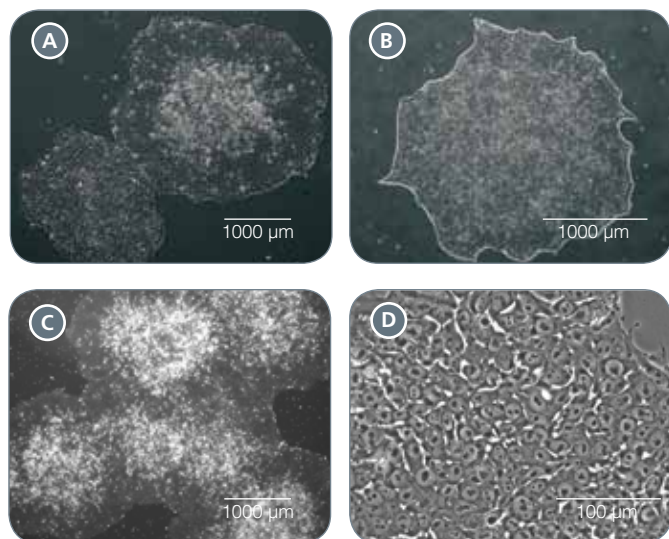


THE CELL EXPERTS™ | [WWW.STEMCELL.COM](http://WWW.STEMCELL.COM)

CATALOG #29064 | VERSION 1.0.0 | DECEMBER 2010

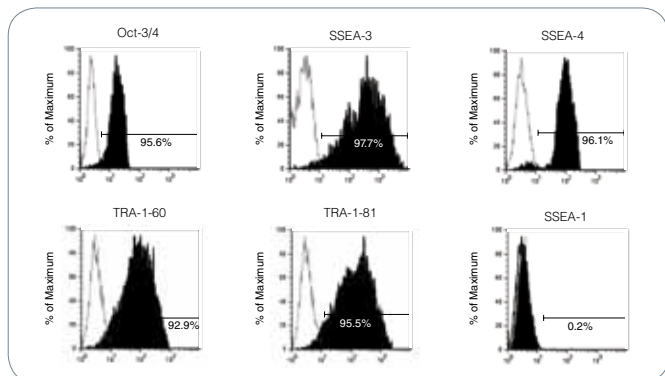
TOLL-FREE T. 1 800 667 0322 • T. +1 604 877 0713 • TOLL-FREE F. 1 800 567 2899 • F. +1 604 877 0704  
[ORDERS@STEMCELL.COM](mailto:ORDERS@STEMCELL.COM) • [INFO@STEMCELL.COM](mailto:INFO@STEMCELL.COM) • FOR FULL CONTACT DETAILS WORLDWIDE VISIT OUR WEBSITE

FOR RESEARCH USE ONLY. NOT FOR THERAPEUTIC OR DIAGNOSTIC USE.

**FIGURE 2.** Morphology of hESCs & hiPSCs Cultured in mTeSR™1


Morphology of hESC and hiPSC lines cultured in mTeSR™1 varies slightly compared to feeder-containing or conditioned medium cultures - colonies have defined edges and high nucleus to cytoplasm ratio. Representative hiPSC lines **(A)** MSC-iPSC1 and **(B)** iPSC(IMR90)-3 and **(C and D)** H9 hESCs are pictured.

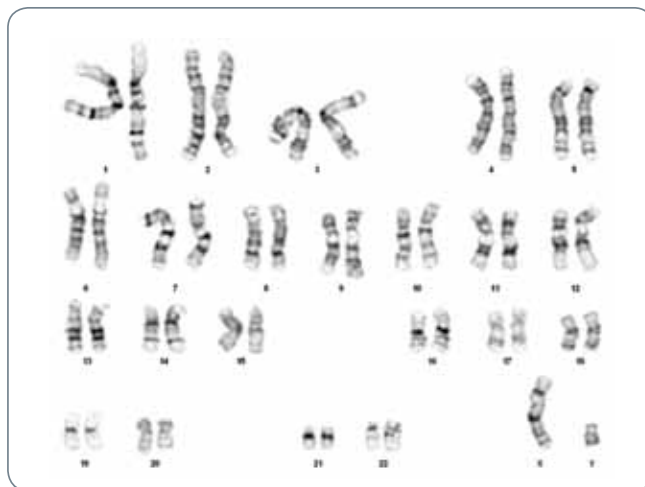
hiPSC photographs courtesy of M. O'Connor and C. Eaves, The Vancouver Human Embryonic Stem Cell Core Facility.

**FIGURE 4.** hESCs & hiPSCs Cultured in mTeSR™1 Retain Expression of Pluripotency Markers


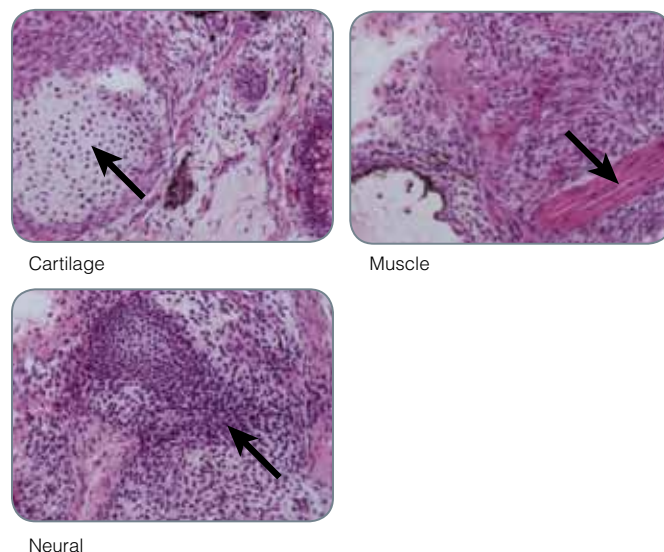
Flow cytometric analysis of H9 hESCs maintained in mTeSR™1 for 17 passages.

## References

- Yu J, Vodyanik MA, Smuga-Otto K, Antosiewicz-Bourget J, Frane JL, Tian S, Nie J, Jonsdottir GA, Ruotti V, Stewart R, Slukvin II, Thomson JA. Induced pluripotent stem cell lines derived from human somatic cells. *Science* 318:1917-1920, 2007
- Ludwig TE, Bergendahl V, Levenstein ME, Yu J, Probasco MD, Thomson JA. Feeder-independent culture of human embryonic stem cells. *Nat Methods* 3:637-646, 2006

**FIGURE 3.** hESCs & hiPSCs Cultured Long-Term in mTeSR™1 Retain Normal Karyotype


Chromosomal analysis of H1 hESCs cultured in mTeSR™1 for 48 passages shows that normal karyotype is retained during long-term passaging.

**FIGURE 5.** hESCs & hiPSCs Cultured in mTeSR™1 are Pluripotent


H9 hESCs were cultured for 6 passages in mTeSR™1 then injected subcutaneously into immunocompromised mice. The resulting teratoma contained cell types from all 3 germ layers. Representative tissue types are shown.