

# MethoCult™ GF M3434



Scientists Helping Scientists™ | WWW.STEMCELL.COM

TOLL FREE PHONE 1 800 667 0322 • PHONE +1 604 877 0713

INFO@STEMCELL.COM • TECHSUPPORT@STEMCELL.COM

FOR GLOBAL CONTACT DETAILS VISIT OUR WEBSITE

## Methylcellulose-based medium with recombinant cytokines (including EPO) for mouse cells

Catalog # 03444      24 x 3 mL  
03434      100 mL

## Product Description

### Complete Methylcellulose-Based Medium for Colony-Forming Unit (CFU) Assays for Mouse Cells

MethoCult™ GF M3434 is optimized for the detection and quantification of mouse hematopoietic progenitor cells in bone marrow, spleen, peripheral blood, and fetal liver samples using CFU assays.

MethoCult™ GF M3434 is formulated to support optimal growth of erythroid progenitor cells (CFU-E and BFU-E), granulocyte-macrophage progenitor cells (CFU-GM, CFU-G, and CFU-M), and multipotential granulocyte, erythroid, macrophage, megakaryocyte progenitor cells (CFU-GEMM).

## Properties

**Storage:** Store at -20°C.

NOTE: Product may be shipped with dry ice or ice packs and may be received thawed.

**Shelf Life:** Stable until expiry date (EXP) on label.

- Contains:**
- Methylcellulose in Iscove's MDM
  - Fetal bovine serum
  - Bovine serum albumin
  - Recombinant human insulin
  - Human transferrin (iron-saturated)
  - 2-Mercaptoethanol
  - Recombinant mouse stem cell factor (SCF)
  - Recombinant mouse interleukin 3 (IL-3)
  - Recombinant human interleukin 6 (IL-6)
  - Recombinant human erythropoietin (EPO)
  - Supplements

This product contains material derived from human plasma. Donors have been tested and found negative for HIV-1 and -2, hepatitis B, and hepatitis C prior to donation. However, this product should be considered potentially infectious and treated in accordance with universal handling precautions.

## Handling/Directions for Use

NOTE: If product is received partially thawed, place immediately at -20°C or thaw and aliquot as described below.

NOTE: Do not use pipettes to dispense methylcellulose as the volume dispensed will not be accurate. Syringes and large bore blunt-end needles should be used for accurate dispensing of viscous methylcellulose medium and to prevent needle-stick injuries.

### A. TO PREPARE 100 ML BOTTLE (Catalog #03434)

1. Thaw MethoCult™ GF M3434 at room temperature (15 - 25°C) or overnight at 2 - 8°C. Do not thaw at 37°C.
2. Shake vigorously for 1 - 2 minutes and then let stand for at least 5 minutes to allow bubbles to rise to the top before aliquoting.
3. Using a luer lock syringe (3 mL [Catalog #28230] or 6 mL) attached to a 16 Gauge Blunt-End Needle (Catalog #28110), aliquot as follows:
  - 3 mL per tube for 1.1 mL duplicate cultures
  - 4 mL per tube for 1.1 mL triplicate cultures

NOTE: Tubes can be used immediately or stored at -20°C. Do not exceed the shelf life of the medium. After thawing aliquoted tubes, mix well and use immediately. Do not re-freeze.

**B. TO PREPARE 3 ML TUBES (Catalog #03444)**

Thaw 3 mL tubes of MethoCult™ GF M3434 at room temperature (15 - 25°C) or overnight at 2 - 8°C. Do not thaw at 37°C. Mix thoroughly.

**NOTE:** After thawing tubes of MethoCult™, use immediately or store at 2 - 8°C for up to 1 month. Do not re-freeze.

For recommended cell plating concentrations, setup of mouse CFU assays, and counting and classification of mouse colonies, refer to the Technical Manual: Mouse Colony-Forming Unit (CFU) Assays Using MethoCult™ (Document #10000005597), available at [www.stemcell.com](http://www.stemcell.com), or contact us to request a copy.

## Related Products

For related products, including specialized culture and storage media, supplements, antibodies, cytokines, and small molecules, visit [www.stemcell.com/HSPCworkflow](http://www.stemcell.com/HSPCworkflow), or contact us at [techsupport@stemcell.com](mailto:techsupport@stemcell.com). For available fresh and cryopreserved peripheral blood, cord blood, and bone marrow products, visit [www.stemcell.com/primarycells](http://www.stemcell.com/primarycells).

## References

Miller CL & Lai B. (2005) Human and mouse hematopoietic colony-forming cell assays. In: Helgason CD & Miller CL (Eds.). Basic Cell Culture Protocols (pp. 71–89). Totowa, New Jersey: Humana Press Inc.

PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED. FOR ADDITIONAL INFORMATION ON QUALITY AT STEMCELL, REFER TO [WWW.STEMCELL.COM/COMPLIANCE](http://WWW.STEMCELL.COM/COMPLIANCE).

Copyright © 2022 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, Scientists Helping Scientists, and MethoCult are trademarks of STEMCELL Technologies Canada Inc. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.