

# FreSR™-S

## Animal component-free medium for cryopreserving ES and iPS cells as single cells

Catalog # 05859 50 mL



Scientists Helping Scientists™ | [WWW.STEMCELL.COM](http://WWW.STEMCELL.COM)

TOLL FREE PHONE 1 800 667 0322 • PHONE +1 604 877 0713

[INFO@STEMCELL.COM](mailto:INFO@STEMCELL.COM) • [TECHSUPPORT@STEMCELL.COM](mailto:TECHSUPPORT@STEMCELL.COM)

FOR GLOBAL CONTACT DETAILS VISIT OUR WEBSITE

## Product Description

FreSR™-S is a defined, serum-free, and animal component-free medium for the cryopreservation of human pluripotent stem cells. This complete and ready-to-use medium is recommended for human embryonic stem (ES) or induced pluripotent stem (iPS) cells cultured in mTeSR™1 (Catalog #85850), mTeSR™ Plus (Catalog #100-0276), TeSR™-E8™ (Catalog #05990), or TeSR™-AOF (Catalog #100-0401). Cryopreserved ES or iPS cells should be stored at -135°C (liquid nitrogen) or colder.

NOTE: This product is not recommended for freezing cell aggregates.

- Defined, serum-free and animal component-free medium for cryopreserving ES/iPS cells as single cells
- Quickly recover ES/iPS cell colonies after thawing
- Reproducibly high recovery rates
- Optimized for cryopreserving ES/iPS cells cultured in TeSR™ maintenance media
- Preserves ES/iPS cell pluripotency and expansion capacities
- Convenient, ready-to-use format

## Properties

**Storage:** Store at 2 - 8°C.

**Shelf Life:** Stable until expiry date (EXP) on label.

Please refer to the Safety Data Sheet (SDS) for hazard information. This product contains components dissolved in dimethyl sulfoxide (DMSO). DMSO is a strong solvent and skin penetrant, and can transport many substances through the skin. DMSO can also penetrate some protective glove materials including latex and silicone. Extra caution should be utilized when handling this product.

## Handling/Directions for Use

NOTE: For complete instructions on culturing ES and iPS cells, on preparing a single-cell suspension using Gentle Cell Dissociation Reagent (Catalog #100-0485), and on coating plates with Vitronectin XF™ (Catalog #07180) or Corning® Matrigel® (Corning Catalog #354277), refer to the Technical Manual: Maintenance of Human Pluripotent Stem Cells in mTeSR™1 (Document #10000005505), mTeSR™ Plus (Document #10000007757), TeSR™-E8™ (Document #10000005516), or TeSR™-AOF (Document #10000008160). These documents are available at [www.stemcell.com](http://www.stemcell.com), or contact us to request a copy.

### CRYOPRESERVING CELLS

1. Prepare a single-cell suspension of human ES or iPS cells using the desired dissociation reagent.
2. Perform a viable cell count using Trypan Blue (Catalog #07050) and a hemocytometer.
3. Centrifuge cells at 300 x g for 5 minutes to obtain a cell pellet.
4. Carefully remove the supernatant with a pipette, leaving a small amount of medium to ensure the cell pellet is not disturbed. Resuspend the cell pellet by gently flicking the tube.
5. Wipe the outside of the FreSR™-S container with 70% ethanol or isopropanol before opening.
6. Add cold (2 - 8°C) FreSR™-S to obtain a cell suspension of  $1 \times 10^6$  cells/mL and mix thoroughly.
7. Transfer 1 mL of the single-cell suspension into each cryovial.
8. Cryopreserve cells using a standard slow rate-controlled cooling protocol (approximately -1°C/minute) or an isopropanol freezing container and store at liquid nitrogen temperature (-135°C).

NOTE: Long-term storage at -80°C is not recommended.

## THAWING CELLS

1. Have all tubes, warmed (15 - 25°C) desired culture medium (e.g. mTeSR™1, mTeSR™ Plus, TeSR™-E8™, or TeSR™-AOF) and DMEM/F-12 with 15 mM HEPES (Catalog #36254), and coated cultureware ready before starting the protocol to ensure that the thawing procedure is done as quickly as possible.
2. Add Y-27632 (Catalog #72302) to culture medium to reach a final concentration of 10 µM.
3. Wipe the outside of the vial of cells with 70% ethanol or isopropanol.
4. In a biosafety cabinet, twist the cap a quarter-turn to relieve internal pressure and then retighten.
5. Quickly thaw cells in a 37°C water bath by gently shaking the vial. Remove the vial when a small frozen cell pellet remains. Do not vortex cells.
6. Wipe the outside of the vial with 70% ethanol or isopropanol.
7. Use a 1 mL pipettor to slowly transfer the contents of the cryovial to a 15 mL conical tube (e.g. Catalog #38009) containing 5 - 7 mL of DMEM/F-12 with 15 mM HEPES.
8. Centrifuge cells at 300 x *g* for 5 minutes at room temperature (15 - 25°C).
9. Carefully remove the supernatant with a pipette, leaving a small amount of medium to ensure the cell pellet is not disturbed.
10. Add 1 mL of culture medium containing 10 µM Y-27632 to the tube. Mix gently.
11. Plate cells onto coated cultureware.  
NOTE: In general, one frozen cryovial containing 1 x 10<sup>6</sup> cells can be thawed and plated into 1 - 2 wells of a 6-well plate.
12. Place the cultureware in a 37°C incubator. Move the cultureware in several quick, short, back-and-forth and side-to-side motions to evenly distribute the cells across the surface.
13. Perform daily medium changes using culture medium without Y-27632, and visually assess cultures to monitor growth until the next passaging time (i.e. 80 - 90% confluent). This takes approximately 2 - 5 days after thawing.  
NOTE: This time may vary when using different cell lines. Cultures should be monitored under the microscope until the optimal passaging time is determined.
14. Passage cultures using standard techniques to generate cell aggregates (e.g. using ReLeSR™ [Catalog #05872]).  
NOTE: It is generally not recommended to perform serial single-cell passaging due to the increased risk of karyotype abnormalities (Buzzard JJ et al.; Draper JS et al.).

## References

1. Buzzard JJ et al. (2004) Karyotype of human ES cells during extended culture. *Nat Biotechnol* 22(4): 381–2.
2. Draper JS et al. (2004) Recurrent gain of chromosomes 17q and 12 in cultured human embryonic stem cells. *Nat Biotechnol* 22(1): 53–4.

PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED. FOR ADDITIONAL INFORMATION ON QUALITY AT STEMCELL, REFER TO [WWW.STEMCELL.COM/COMPLIANCE](http://WWW.STEMCELL.COM/COMPLIANCE).

Copyright © 2023 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, Scientists Helping Scientists, ReLeSR, and FreSR are trademarks of STEMCELL Technologies Canada Inc. E8, mTeSR, and TeSR are trademarks of WARF. Vitronectin XF is developed and manufactured by Nucleus Biologics, and Vitronectin XF is a trademark of Nucleus Biologics. Corning and Matrigel are registered trademarks of Corning Incorporated. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.