

# DNase I Solution (1 mg/mL)

## Cell dissociation reagent

Catalog # 07900                      1 mL  
100-0762                              10 x 1 mL



Scientists Helping Scientists™ | [WWW.STEMCELL.COM](http://WWW.STEMCELL.COM)

TOLL FREE PHONE 1 800 667 0322 • PHONE +1 604 877 0713

[INFO@STEMCELL.COM](mailto:INFO@STEMCELL.COM) • [TECHSUPPORT@STEMCELL.COM](mailto:TECHSUPPORT@STEMCELL.COM)

FOR GLOBAL CONTACT DETAILS VISIT OUR WEBSITE

## Product Description

Deoxyribonuclease (DNase) I Solution (1 mg/mL) is useful for reducing or preventing aggregation of concentrated and/or cryopreserved cell suspensions following thawing. The activity range is  $\geq 2000$  units per mg of protein.

## Properties

- Storage:** Store at -20°C.
- Shelf Life:** Stable until expiry date (EXP) on label.
- Contains:**
- Bovine pancreatic deoxyribonuclease I (1 mg/mL)
  - Phosphate-buffered saline

## Handling/Directions for Use

Storage at 2 - 8°C is not recommended. Use the protocol below when cell samples appear to be aggregating; this can occur when cells have been frozen and thawed or after enzymatic tissue digestion. For optimal performance, use cells diluted in a medium that does not contain a chelator of calcium and magnesium ions (e.g. EDTA).

1. Thaw DNase I Solution at room temperature (15 - 25°C) or overnight at 2 - 8°C.
2. Centrifuge cells and carefully remove the supernatant.
3. Resuspend cells in 0.1 mg/mL of DNase I Solution.
4. Incubate at room temperature for 15 minutes.

NOTE: For optimal cell separation results, filter aggregated suspensions through a 37  $\mu$ m Reversible Strainer (Catalog #27215/27250), then resuspend at the appropriate cell concentration in desired medium.

NOTE: For downstream applications that are sensitive to DNase, wash cells with appropriate DNase-free medium.

## Notes

### ACTIVITY UNITS

1 unit refers to the amount of DNase I required to act on 1 mg/mL of DNA (pH 5.0, 25°C) to produce an increase in absorbance of 0.001 per minute at a wavelength of 260 nm.

**This product is hazardous. Please refer to the Safety Data Sheet (SDS).**

PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED.

Copyright © 2022 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, and Scientists Helping Scientists are trademarks of STEMCELL Technologies Canada Inc. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.