

Gentle Collagenase/Hyaluronidase



Scientists Helping Scientists™ | WWW.STEMCELL.COM

TOLL FREE PHONE 1 800 667 0322 • PHONE +1 604 877 0713

INFO@STEMCELL.COM • TECHSUPPORT@STEMCELL.COM

FOR GLOBAL CONTACT DETAILS VISIT OUR WEBSITE

10X Gentle collagenase/hyaluronidase in DMEM

Catalog # 07919

10 mL

Product Description

10X Collagenase/Hyaluronidase in Dulbecco's Modified Eagle's Medium (DMEM) for the gentle enzymatic dissociation of mouse mammary tissue.

Properties

- Storage:** Store at -20°C.
- Shelf Life:** Stable until expiry date (EXP) on label.
- Contains:**
- 10 mg/mL Collagenase
 - 1000 U/mL Hyaluronidase
 - DMEM (1000 mg D-glucose/L)

Please refer to the Safety Data Sheet (SDS) for hazard information.

Materials Required But Not Included

- Ammonium Chloride Solution (Catalog #07800)
- EpiCult™-B Mouse Medium Kit (Catalog #05610)
- Fetal bovine serum (FBS)
- HBSS with 10 mM HEPES, Without Phenol Red (Catalog #37150)
- Gentamicin

Handling / Directions For Use

NOTE: Avoid the use of glass pipettes and tubes when handling mammary epithelial cells as these cells will stick to glass.

DISSOCIATION OF MOUSE MAMMARY TISSUE

1. Thaw Gentle Collagenase/Hyaluronidase at room temperature (15 - 25°C) or overnight at 2 - 8°C.
NOTE: Once thawed, use immediately or aliquot and store at -20°C until the expiry date as indicated on the label. After thawing the aliquots, use immediately. Do not re-freeze.
2. Dilute Gentle Collagenase/Hyaluronidase as outlined below for either 2-hour or overnight dissociation.
NOTE: Approximately 2 - 5 mL of the EpiCult™-B Medium/Collagenase/Hyaluronidase/FBS/gentamicin solution will be required for every 2 mammary glands to be dissociated. Alternatively, mammary glands can be dissociated in DMEM/F-12 with 15 mM HEPES (Catalog #36254) supplemented with 50 µg/mL gentamicin to avoid influences of exogenous growth factors and FBS; however, this may result in lower total viable cell yields.

2-HOUR DISSOCIATION

Dilute 1 part Gentle Collagenase/Hyaluronidase with 4 parts complete EpiCult™-B Medium (Mouse) supplemented with 5% FBS and 50 µg/mL gentamicin in a 15 mL or 50 mL tube (e.g. Catalog #38009 or 38010). Continue to step 3.

OVERNIGHT DISSOCIATION

Dilute 1 part Gentle Collagenase/Hyaluronidase with 9 parts complete EpiCult™-B Medium (Mouse) supplemented with 5% FBS and 50 µg/mL gentamicin in a 15 mL or 50 mL tube (e.g. Catalog #38009 or 38010). Continue to step 4.

3. For 2-hour dissociation, resect mammary glands and transfer to a sterile glass Petri dish. Mince with scalpels in a crosswise pattern until glands are rendered to a paste. This step is not necessary for overnight dissociation.

NOTE: It is essential that the glands are well minced, or total viable epithelial cell yield will be low.

4. Transfer the mammary tissue to the tube containing EpiCult™-B Medium /Collagenase/Hyaluronidase/FBS/gentamicin.

2-HOUR DISSOCIATION: Incubate at 37°C on a rotary/rocking shaker set at 90 rpm for 2 hours. After 1 hour and at the end of incubation, pipette the suspension up and down 20 times using a 1 mL pipettor.

OVERNIGHT DISSOCIATION: Incubate at 37°C for 15 hours (overnight). It is not necessary to use a shaker.

5. Centrifuge the cells at 350 x *g* for 5 minutes and discard the supernatant.
6. Resuspend the cell pellet in a mixture of 1 part cold HBSS with 10 mM HEPES, Without Phenol Red supplemented with 2% FBS and 4 parts Ammonium Chloride Solution and centrifuge at 350 x *g* for 5 minutes. The resultant pellet contains epithelial cell organoids as well as stromal cells and lymphocytes.

To generate a single-cell suspension of mammary epithelial cells, refer to the Technical Bulletin: Monolayer Culture of Mouse Mammary Epithelial Cells (Document #29179), available at www.stemcell.com or contact us to request a copy.

STEMCELL TECHNOLOGIES INC.'S QUALITY MANAGEMENT SYSTEM IS CERTIFIED TO ISO 13485. PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED.

Copyright © 2019 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, Scientists Helping Scientists, and EpiCult are trademarks of STEMCELL Technologies Canada Inc. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.