

## EasySep™ Human TCR Alpha/Beta Depletion Kit

Negative Selection  
Catalog #17847

For processing 1 x 10<sup>9</sup> cells



Scientists Helping Scientists™ | [WWW.STEMCELL.COM](http://WWW.STEMCELL.COM)

TOLL FREE PHONE 1 800 667 0322 • PHONE +1 604 877 0713  
INFO@STEMCELL.COM • TECHSUPPORT@STEMCELL.COM  
FOR GLOBAL CONTACT DETAILS VISIT OUR WEBSITE

Document #1000005219 | Version 00

## Description

Deplete human T cell receptor alpha/beta+ (TCR $\alpha\beta$ +) cells from leukapheresis samples or expanded TCR $\alpha\beta$  knockout cell cultures.

- Fast and easy-to-use
- No columns required
- Isolated cells are untouched

This kit targets TCR $\alpha\beta$  cells for removal with an antibody recognizing the TCR $\alpha\beta$  surface marker. Unwanted cells are labeled with antibody and magnetic particles, and separated without columns using an EasySep™ magnet. Desired cells are simply poured off into a new tube. Isolated cells are immediately available for downstream applications such as flow cytometry, culture, or cryopreservation.

## Component Descriptions

COMPONENT NAME	COMPONENT #	QUANTITY	STORAGE	SHELF LIFE	FORMAT
EasySep™ Human TCR Alpha/Beta Depletion Cocktail	17847C	1 x 1 mL	Store at 2 - 8°C. Do not freeze.	Stable until expiry date (EXP) on label.	A combination of monoclonal antibodies in PBS.
EasySep™ Dextran RapidSpheres™ 50103	50103	2 x 1 mL	Store at 2 - 8°C. Do not freeze.	Stable until expiry date (EXP) on label.	A suspension of magnetic particles in water.

PBS - phosphate-buffered saline

Components may be shipped at room temperature (15 - 25°C) but should be stored as indicated above.

## Sample Preparation

For generating a TCR $\alpha\beta$  knockout cell population, refer to the Technical Bulletin: Genome Editing of Human Primary T Cells (Document #27155), available at [www.stemcell.com](http://www.stemcell.com) or contact us to request a copy.

### EXPANDED TCR $\alpha\beta$ KNOCKOUT CELLS

Harvest cultured cells and remove the culture medium by centrifugation. Resuspend the cells at 5 x 10<sup>7</sup> cells/mL in recommended medium.

### LEUKAPHERESIS

1. Add 3 parts Ammonium Chloride (Catalog #07800) to 1 part leukapheresis sample.
2. Incubate on ice for 15 minutes.
3. Centrifuge at 300 x g for 10 minutes at room temperature (15 - 25°C). Carefully remove the supernatant.
4. Wash the cells by topping up the tube with recommended medium. Centrifuge cells at 150 x g for 10 minutes at room temperature with the brake off. Carefully remove the supernatant.
5. Resuspend cells at 5 x 10<sup>7</sup> cells/mL in recommended medium.



## Recommended Medium

EasySep™ Buffer (Catalog #20144), or PBS containing 2% fetal bovine serum (FBS) and 1 mM EDTA. Medium should be free of Ca<sup>++</sup> and Mg<sup>++</sup>.

**Directions for Use – Manual EasySep™ Protocols**

See page 1 for Sample Preparation and Recommended Medium. Refer to Tables 1 and 2 for detailed instructions regarding the EasySep™ procedure for each magnet.


**Table 1. EasySep™ Human TCR Alpha/Beta Depletion Kit Protocol**

		EASYSEP™ MAGNETS	
STEP	INSTRUCTIONS	 <b>EasySep™</b> (Catalog #18000)	<b>“The Big Easy”</b> (Catalog #18001) 
1	Prepare sample at the indicated cell concentration within the volume range.	5 x 10 <sup>7</sup> cells/mL 0.25 - 2 mL	5 x 10 <sup>7</sup> cells/mL 0.5 - 8.5 mL
	Add sample to required tube.	5 mL (12 x 75 mm) polystyrene round-bottom tube (e.g. Catalog #38007)	14 mL (17 x 100 mm) polystyrene round-bottom tube (e.g. Catalog #38008)
2	Add Depletion Cocktail to sample.	50 µL/mL of sample	50 µL/mL of sample
	Mix and incubate.	RT for 5 minutes	RT for 5 minutes
3	Vortex RapidSpheres™. NOTE: Particles should appear evenly dispersed.	30 seconds	30 seconds
4	Add RapidSpheres™ to sample.	50 µL/mL of sample	50 µL/mL of sample
	Mix and incubate.	RT for 3 minutes	RT for 3 minutes
5	Add recommended medium to top up the sample to the indicated volume. Mix by gently pipetting up and down 2 - 3 times.	Top up to 2.5 mL	<ul style="list-style-type: none"> <li>• Top up to 5 mL for samples ≤ 4 mL</li> <li>• Top up to 10 mL for samples &gt; 4 mL</li> </ul>
	Place the tube (without lid) into the magnet and incubate.	RT for 5 minutes	RT for 5 minutes
6	Pick up the magnet, and in one continuous motion invert the magnet and tube,* pouring the depleted cell suspension into a new tube.	Isolated cells are ready for use	Isolated cells are ready for use

RT - room temperature (15 - 25°C)

\* Leave the magnet and tube inverted for 2 - 3 seconds, then return upright. Do not shake or blot off any drops that may remain hanging from the mouth of the tube.

**Table 2. EasySep™ Human TCR Alpha/Beta Depletion Kit Protocol**

		EASYSEP™ MAGNETS	
STEP	INSTRUCTIONS	Easy 50 (Catalog #18002)	
1	Prepare sample at the indicated cell concentration within the volume range.	5 x 10 <sup>7</sup> cells/mL 5 - 40 mL	
	Add sample to required tube.	50 mL (30 x 115 mm) conical tube (e.g. Catalog #38010)	
2	Add Depletion Cocktail to sample.	50 µL/mL of sample	
	Mix and incubate.	RT for 10 minutes	
3	Vortex RapidSpheres™. NOTE: Particles should appear evenly dispersed.	30 seconds	
4	Add RapidSpheres™ to sample.	100 µL/mL of sample	
	Mix and incubate.	RT for 5 minutes	
5	Add recommended medium to top up sample to the indicated volume. Mix by gently pipetting up and down 2 - 3 times.	<ul style="list-style-type: none"> <li>• Top up to 25 mL for samples ≤ 20 mL</li> <li>• Top up to 50 mL for samples &gt; 20 mL</li> </ul>	
	Place the tube (without lid) into the magnet and incubate.	RT for 10 minutes	
6	Carefully pipette (do not pour) the depleted cell suspension into a new tube.	Use a new 50 mL tube	
7	Remove the tube from the magnet and place the new tube (without lid) into the magnet and incubate for a second separation.	RT for 10 minutes (total of 2 x 10-minute separations)	
8	Carefully pipette (do not pour) the depleted cell suspension into a new tube.	Isolated cells are ready for use	

RT - room temperature (15 - 25°C)

## Notes and Tips


### ASSESSING PURITY

For purity assessment of residual TCRαβ+ cells by flow cytometry, use the following fluorochrome-conjugated antibody clones:

- Anti-human TCRαβ antibody, clone IP26 (partially blocked), or
- Anti-human TCRαβ antibody, clone T10B9.1A31 (partially blocked)

NOTE: The expression of endogenous TCRαβ on human primary T cells can be disrupted using ArciTect™ CRISPR/Cas9-mediated gene modification. The TCRαβ-knockout efficiency typically ranges from 60 - 90%. The expression level of TCRαβ in the starting population from expanded TCRαβ knockout cell culture varies with the TCRαβ knockout efficiency.

## Data

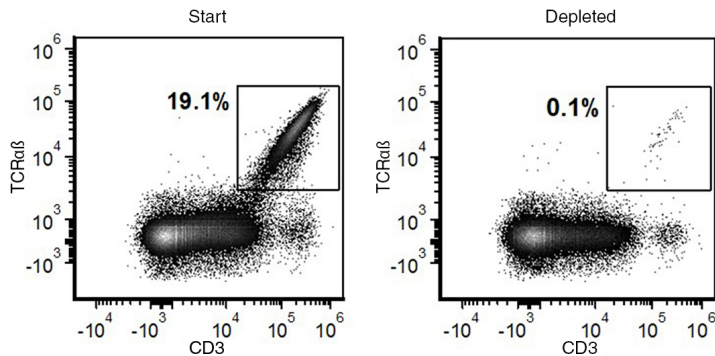

In the above example, the frequencies of CD3+TCRαβ+ cells in the starting and depleted fractions are 19.1% and 0.1%, respectively.

STEMCELL TECHNOLOGIES INC.'S QUALITY MANAGEMENT SYSTEM IS CERTIFIED TO ISO 13485. PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED.

Copyright © 2019 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, Scientists Helping Scientists, ArciTect, EasySep, and RapidSpheres are trademarks of STEMCELL Technologies Inc. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.