

Conditional Reprogramming (CR) Medium



Scientists Helping Scientists™ | WWW.STEMCELL.COM

TOLL FREE PHONE 1 800 667 0322 • PHONE +1 604 877 0713

INFO@STEMCELL.COM • TECHSUPPORT@STEMCELL.COM

FOR GLOBAL CONTACT DETAILS VISIT OUR WEBSITE

Catalog #100-0352

250 mL

Conditional reprogramming medium for epithelial cell expansion

Product Description

Conditional Reprogramming (CR) Medium is optimized to use with 3T3-J2 Irradiated Feeder Cells (Catalog #100-0353) to expand epithelial cells without accumulating karyotypic modifications. Epithelial cells seeded into CR Medium revert to a stem-like state that promotes long-term expansion, while maintaining differentiation potential (Liu X et al. 2012; Liu X et al. 2017). This conditional proliferative capability is deactivated once the medium is removed from the cells (Liu X et al. 2012; Liu X et al. 2017). CR Medium is suitable to propagate primary epithelial cells from healthy or tumor tissues (airway, retina, prostate, breast, intestine, pancreas, liver biliary duct, etc.; Liu X et al. 2012; Liu X et al. 2017; Suprynowicz FA et al.) from multiple species, including humans, mouse and rat. This product is free of antibiotics and intended for Research Use Only.

NOTE: CR Medium must be supplemented with cholera toxin prior to use.

Properties

Storage: Store at -20°C.

Shelf Life: Stable for 2 years from date of manufacture (MFG) on label. Protect from prolonged exposure to light.

Materials Required But Not Included

PRODUCT NAME	CATALOG #
Cholera toxin	e.g. Sigma C8052
3T3-J2 Irradiated Feeder Cells	100-0353

Preparation of Complete CR Medium

Use sterile technique to prepare complete CR Medium (CR Medium + cholera toxin). The following example is for preparing 250 mL of complete CR Medium. If preparing other volumes, adjust accordingly.

1. Thaw CR Medium at room temperature (15 - 25°C) or overnight at 2 - 8°C. Mix thoroughly.

NOTE: Once thawed, use immediately or store at 2 - 8°C for up to 1 month.

2. Prepare cholera toxin stock solution by reconstituting cholera toxin in CR Medium to a final concentration of 1 mg/mL. Mix thoroughly.

NOTE: If not used immediately, aliquot and store at 2 - 8°C for up to 6 months.

3. Add the appropriate volume of cholera toxin stock solution (1 mg/mL; prepared in step 2) to 250 mL of CR Medium to achieve a final concentration of 8.6 - 20 ng/mL.

For example, add 2.2 µL of cholera toxin stock solution (1 mg/mL) for a final concentration of 8.6 ng/mL.

Mix thoroughly.

NOTE: Complete CR Medium does not contain antibiotics; add antibiotics if desired.

NOTE: If not used immediately, store complete CR Medium at 2 - 8°C for up to 1 month. Do not exceed the shelf life of the individual components.

Directions For Use

CO-CULTURE OF EPITHELIAL CELLS WITH 3T3-J2 IRRADIATED FEEDER CELLS

The following protocol is for expanding cultured primary epithelial cells in a single T-25 cm² flask using 3T3-J2 Irradiated Feeder Cells (Catalog #100-0353) and complete CR Medium. If using other cultureware, adjust cell numbers and volumes accordingly.

NOTE: In this protocol, 3T3-J2 Irradiated Feeder Cells and epithelial cells are seeded at the same time.

1. Prepare complete CR Medium (see Preparation section). Warm in a 37°C water bath.

2. Thaw and prepare 3T3-J2 Irradiated Feeder Cells. For complete instructions, refer to the applicable Product Information Sheet (Document #10000007812), available at www.stemcell.com or contact us to request a copy.
3. Add 7.5×10^5 freshly thawed and washed 3T3-J2 Irradiated Feeder Cells to 5 mL of complete CR Medium. This number of feeder cells will result in a final seeding density of 3×10^4 viable cells/cm² (a range of $2.5 - 4 \times 10^4$ viable cells/cm² is recommended).
4. Add the desired number of epithelial cells to the feeder cell suspension.
5. Add the cell suspension to the T-25 cm² flask and incubate at 37°C and 5% CO₂. Perform a full-medium change every 2 - 3 days.

Related Products

For a complete list of related products available from STEMCELL Technologies, including specialized cell culture and storage media, antibodies, cytokines, and small molecules, visit www.stemcell.com or contact us at techsupport@stemcell.com.

References

- Liu X et al. (2017) Conditional reprogramming and long-term expansion of normal and tumor cells from human biospecimens. *Nat Protoc* 12(2): 439–51.
- Liu X et al. (2012) ROCK inhibitor and feeder cells induce the conditional reprogramming of epithelial cells. *Am J Pathol* 180(2): 599–607.
- Suprynowicz FA et al. (2012) Conditionally reprogrammed cells represent a stem-like state of adult epithelial cells. *Proc Natl Acad Sci USA* 109(49): 20035–40.

Conditional Reprogramming (CR) Medium is manufactured by Propagenix Inc.

PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED.

Copyright © 2020 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, and Scientists Helping Scientists are trademarks of STEMCELL Technologies Canada Inc. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.