




4721 Calle Carga  
Camarillo, CA 93012  
[www.pbsbiotech.com](http://www.pbsbiotech.com)

VERTICAL-WHEEL®  
BIOREACTORS



**PBS**MINI

PBS 0.1 and 0.5 Single-Use Vessels

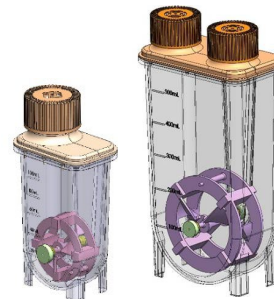

**Quick Guide**

IC01024 Rev A

## Highlights for PBS MINI

The PBS 0.1 and PBS 0.5 Single-Use Vessels are primarily constructed from injection-molded polycarbonate and are sterilized by gamma irradiation to be ready-for-use by the customer.

1. The MINI Base can be used with either PBS 0.1 or PBS 0.5 vessels.
2. Working volume range can be found in the MINI Base User Manual or on our website [www.pbsbiotech.com](http://www.pbsbiotech.com).
3. Do not press the buttons on the MINI Base when the display is ON.



## PBS MINI Vessel Handling

1. Do not re-use vessel. Vessels are designed to be single-use disposables.
2. Do not overtighten the vent cap(s).
3. Ensure the cap(s) are gently hand-tightened between uses to minimize contamination risks.
4. Do not compress the vessel from the sides.