

# Agar Leukocyte Conditioned Medium



**Source of colony-stimulating factors for assays of human hematopoietic progenitor cells**

Catalog # 02300      25 mL

Scientists Helping Scientists™ | [WWW.STEMCELL.COM](http://WWW.STEMCELL.COM)

TOLL FREE PHONE 1 800 667 0322 • PHONE +1 604 877 0713

[INFO@STEMCELL.COM](mailto:INFO@STEMCELL.COM) • [TECHSUPPORT@STEMCELL.COM](mailto:TECHSUPPORT@STEMCELL.COM)

FOR GLOBAL CONTACT DETAILS VISIT OUR WEBSITE

## Product Description

Agar Leukocyte Conditioned Medium (Agar-LCM) is a source of colony-stimulating factors for assays of human hematopoietic progenitor cells from light density fractions of bone marrow and peripheral blood. It is suitable for the growth of colony-forming unit erythroid (CFU-E), burst-forming unit-erythroid (BFU-E), colony-forming unit-macrophage (CFU-M), colony-forming unit-granulocyte (CFU-G), colony-forming unit-granulocyte, macrophage (CFU-GM) and colony-forming unit-granulocyte, erythroid, macrophage, megakaryocyte (CFU-GEMM) when added to MethoCult™ medium at a final concentration of 10% with the addition of erythropoietin (EPO; Catalog #02625). Agar-LCM is prepared using normal human peripheral blood leukocytes in an agar-containing medium.

## Properties

- Storage:** Store at -20°C.
- Shelf Life:** Stable until expiry date (EXP) on label.
- Contains:**
- Agar Leukocyte Conditioned Medium
  - Fetal bovine serum (10%)
  - Iscove's MDM
  - $1 \times 10^4$  2-Mercaptoethanol (ME)

Please refer to the Safety Data Sheet (SDS) for hazard information.

## Handling / Directions For Use

Thaw Agar-LCM at room temperature (15 - 25°C) or overnight at 2 - 8°C.

NOTE: After thawing, some particulate matter may appear; the product may be clarified by centrifugation. There should be no loss in bioactivity as a result of this procedure. Keep on ice while in use.

Preparation of MethoCult™ Methylcellulose-based Medium Containing Agar-LCM for Detection of Human Hematopoietic Progenitor Cells (BFU-E, CFU-GM, CFU-GEMM):

1. Thaw 80 mL bottle of MethoCult™ H4230 (Catalog #04230) at room temperature (15 - 25°C) or overnight at 2 - 8°C. Mix contents by vigorous shaking.
2. Add 10 mL of Agar-LCM. Add EPO to yield a final concentration of 3 U/mL. Add Iscove's Modified Dulbecco's Medium (IMDM; Catalog #36150) to give a total volume of 100 mL.
3. Mix medium by vigorous shaking for 1 - 2 minutes. Let stand for 5 minutes to allow bubbles to rise to the top.
4. Dispense 3 mL of medium into sterile culture tubes using a 6 or 12 mL syringe attached to a 16 Gauge Blunt-End Needle (Catalog #28110).

NOTE: If not used immediately, cap tubes tightly and store at -20°C. Thaw tubes at room temperature (15 - 25°C).

5. Prepare a suspension of human cells in Iscove's MDM with 2% FBS (Catalog #07700) at 10X the final concentration required for plating.
6. Add 0.3 mL of cells to 3.0 mL of MethoCult™ medium.  
NOTE: The addition of cells at a 1:10 (v/v) ratio to methylcellulose-based medium maintains optimal medium viscosity.
7. Vortex tubes to ensure contents are thoroughly mixed. Let stand for 2 - 3 minutes to allow bubbles to rise to the top.
8. Dispense 1.1 mL of MethoCult™ mixture containing cells into each of two 35 mm Culture Dishes (Catalog #27100) using a 3 mL syringe attached to a 16 Gauge Blunt-End Needle.
9. Incubate for 12 - 14 days at 37°C in 5% CO<sub>2</sub> with ≥ 95% humidity.
10. Score colonies using an inverted microscope.

NOTE: For additional information refer to the Technical Manual: Human Colony-Forming Unit (CFU) Assays Using MethoCult™ (Document #28404), available on our website at [www.stemcell.com](http://www.stemcell.com) or contact us to request a copy.

STEMCELL TECHNOLOGIES INC.'S QUALITY MANAGEMENT SYSTEM IS CERTIFIED TO ISO 13485. PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED.

Copyright © 2016 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, MethoCult, and Scientists Helping Scientists are trademarks of STEMCELL Technologies Canada Inc. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.