

B. TO PREPARE 3 ML TUBES (Catalog # 03444)

Thaw 3 mL tubes of MethoCult™ GF M3434 at room temperature (15 - 25°C) or overnight at 2 - 8°C. Do not thaw MethoCult™ at 37°C. Mix thoroughly.

NOTE: After thawing tubes of MethoCult™, use immediately or store at 2 - 8°C for up to 1 month. Do not re-freeze.

For recommended cell plating concentrations, setup of mouse CFU assays, and counting and classification of mouse colonies, refer to the Technical Manual: Mouse Colony-Forming Unit (CFU) Assays Using MethoCult™ (Document #28405), available at www.stemcell.com or contact us to request a copy.

Notes and Tips

RELATED PRODUCTS

For related products, including specialized culture and storage media, supplements, antibodies, cytokines, and small molecules, visit www.stemcell.com/HSPCworkflow or contact us at techsupport@stemcell.com.

References

Miller CL & Lai B. (2005) Human and mouse hematopoietic colony-forming cell assays. In: Helgason CD & Miller CL (Eds.). Basic Cell Culture Protocols (pp. 71–89). Totowa, New Jersey: Humana Press Inc.

STEMCELL TECHNOLOGIES INC.'S QUALITY MANAGEMENT SYSTEM IS CERTIFIED TO ISO 13485. PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED.

Copyright © 2018 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, Scientists Helping Scientists, and MethoCult are trademarks of STEMCELL Technologies Canada Inc. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.