

MethoCult™ M3630

Methylcellulose-based medium for mouse pre-B lymphoid progenitor cells

Catalog # 03630 100 mL



Scientists Helping Scientists™ | WWW.STEMCELL.COM

TOLL FREE PHONE 1 800 667 0322 • PHONE +1 604 877 0713

INFO@STEMCELL.COM • TECHSUPPORT@STEMCELL.COM

FOR GLOBAL CONTACT DETAILS VISIT OUR WEBSITE

Product Description

Complete Methylcellulose-Based Medium for Colony-Forming Unit (CFU) Assays for Mouse Cells

MethoCult™ M3630 is optimized for the detection and quantification of mouse pre-B progenitor cells in bone marrow and Whitlock-Witte long-term cultures using CFU assays.

MethoCult™ M3630 is formulated to support optimal growth of CFU pre-B lymphoid progenitor cells.

Properties

- Storage:** Store at -20°C.
- Shelf Life:** Stable until expiry date (EXP) on label.
- Contains:**
- Methylcellulose in Iscove's MDM
 - Fetal bovine serum
 - 2-Mercaptoethanol
 - Recombinant human interleukin 7 (IL-7)
 - Supplements

Handling / Directions For Use

NOTE: If product is received partially thawed, place immediately at -20°C or thaw and aliquot as described below. Do not use MethoCult™ past the expiry date as indicated on the label.

NOTE: Do not use pipettes to dispense methylcellulose as the volume dispensed will not be accurate. Syringes and large bore blunt-end needles should be used for accurate dispensing of viscous methylcellulose medium and to prevent needle-stick injuries.

1. Thaw 100 mL bottle of MethoCult™ M3630 at room temperature (15 - 25°C) or overnight at 2 - 8°C. Do not thaw methylcellulose-based media at 37°C.
2. Shake vigorously for 1 - 2 minutes and then let stand for at least 5 minutes to allow bubbles to rise to the top before aliquoting.
3. Using a luer lock syringe (3 mL [Catalog #28230] or 6 mL) attached to a 16 gauge Blunt-End Needle (Catalog #28110), aliquot as follows:
 - 3 mL per tube for 1.1 mL duplicate cultures
 - 4 mL per tube for 1.1 mL triplicate cultures

NOTE: Tubes can be used immediately or stored at -20°C. After thawing aliquoted tubes of MethoCult™, mix well and use immediately. Do not re-freeze.

For recommended cell plating concentrations, setup of mouse CFU assays, and counting and classification of mouse colonies, refer to the Technical Manual: Mouse Colony-Forming Unit (CFU) Assays Using MethoCult™ (Document #28405), available at www.stemcell.com or contact us to request a copy.

Notes and Tips

RELATED PRODUCTS

For related products, including specialized culture and storage media, supplements, antibodies, cytokines, and small molecules, visit www.stemcell.com/HSPCworkflow or contact us at techsupport@stemcell.com.

References

Miller CL & Lai B. (2005) Human and mouse hematopoietic colony-forming cell assays. In: Helgason CD & Miller CL (Eds.). Basic Cell Culture Protocols (pp. 71–89). Totowa, New Jersey: Humana Press Inc.

STEMCELL TECHNOLOGIES INC.'S QUALITY MANAGEMENT SYSTEM IS CERTIFIED TO ISO 13485. PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED.

Copyright © 2018 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, Scientists Helping Scientists, and MethoCult are trademarks of STEMCELL Technologies Canada Inc. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.