

# MyeloCult™ H5100

## Medium for long-term culture of human cells

Catalog #	05100	100 mL
	05150	500 mL



Scientists Helping Scientists™ | [WWW.STEMCELL.COM](http://WWW.STEMCELL.COM)

TOLL FREE PHONE 1 800 667 0322 • PHONE +1 604 877 0713

[INFO@STEMCELL.COM](mailto:INFO@STEMCELL.COM) • [TECHSUPPORT@STEMCELL.COM](mailto:TECHSUPPORT@STEMCELL.COM)

FOR GLOBAL CONTACT DETAILS VISIT OUR WEBSITE

## Product Description

MyeloCult™ H5100 is optimized for initiation and maintenance of myeloid long-term cultures of human hematopoietic cells and stromal cell “feeder” layers. The sera used in this formulation have been pre-tested and selected for their ability to support long-term myelopoiesis by primitive human hematopoietic cells (e.g. in long-term culture-initiating cell [LTC-IC] assays).

MyeloCult™ H5100 requires the addition of freshly prepared Hydrocortisone (Catalog #07904).

## Properties

<b>Storage:</b>	Store at -20°C.
<b>Shelf Life:</b>	Stable until expiry date (EXP) on label.
<b>Contains:</b>	<ul style="list-style-type: none"><li>• Horse serum</li><li>• Fetal bovine serum</li><li>• 2-Mercaptoethanol</li><li>• Minimum Essential Medium (MEM) Alpha</li><li>• Supplements</li></ul>

## Handling / Directions For Use

For detailed instructions refer to the Technical Manual: Human Long-Term Culture-Initiating Cell (LTC-IC) Assay (Document #28412), available on our website at [www.stemcell.com](http://www.stemcell.com) or contact us to request a copy.

1. Thaw MyeloCult™ H5100 medium in refrigerator (2 - 8°C), at room temperature (15 - 25°C), or at 37°C until just thawed. Mix well by swirling.  
NOTE: If not used immediately, store at 2 - 8°C for up to 1 month.  
NOTE: If precipitate or crystals are present after thawing, allow the medium to equilibrate to room temperature (15 - 25°C) or place at 37°C for up to 1 hour and mix well by swirling. If precipitate or crystals persist, centrifuge or filter using a 0.2 µm low protein binding filter before use.
2. Add freshly prepared and filter-sterilized hydrocortisone to give a final concentration of 10<sup>-6</sup> M.  
NOTE: After addition of hydrocortisone, the medium is stable for up to 1 week when stored at 2 - 8°C.
3. Cultures should be established in tissue culture-treated plasticware to promote cell attachment and adherent layer formation.  
NOTE: During medium changes, care should be taken to avoid disturbing the adherent layer. Do not touch the bottom of the dish when pipetting; instead, rest the pipette against the side of the dish. When dispensing, allow medium to run slowly down the side of the dish rather than directly onto the surface of the adherent layer.

## Notes and Tips

### RELATED PRODUCTS

For related products, including specialized culture and storage media, supplements, antibodies, cytokines, and small molecules, visit [www.stemcell.com/HSPCworkflow](http://www.stemcell.com/HSPCworkflow) or contact us at [techsupport@stemcell.com](mailto:techsupport@stemcell.com). For available fresh and cryopreserved peripheral blood, cord blood, and bone marrow products in your region, visit [www.stemcell.com/primarycells](http://www.stemcell.com/primarycells).

## References

- Eaves C, Cashman J & Eaves A. (1991) Methodology of long term culture of human hemopoietic cells. *J Tissue Cult Methods* 13: 55–62.
- Hogge DE et al. (1996) Enhanced detection, maintenance, and differentiation of primitive human hematopoietic cells in cultures containing murine fibroblasts engineered to produce human steel factor, interleukin-3, and granulocyte colony-stimulating factor. *Blood* 88(10): 3765–73.
- Miller CL & Eaves CJ. (2002) Long-term culture-initiating cell assays for human and murine cells. *Methods Mol Med* 63: 123–41.
- Petzer AL et al. (1996) Self-renewal of primitive human hematopoietic cells (long-term-culture-initiating cells) in vitro and their expansion in defined medium. *Proc Natl Acad Sci U S A* 93(4): 1470–4.
- Sutherland HJ et al. (1990) Functional characterization of individual human hematopoietic stem cells cultured at limiting dilution on supportive marrow stromal layers. *Proc Natl Acad Sci U S A* 87(9): 3584–8.
- Sutherland HJ, Eaves AC & Eaves CJ. (1991) Quantitative assays for human hemopoietic progenitor cells. In: Gee AP (Ed.), *Bone Marrow Processing and Purging: A Practical Guide*. (pp. 155–71). Boca Raton, FL: CRC Press Inc.

STEMCELL TECHNOLOGIES INC.'S QUALITY MANAGEMENT SYSTEM IS CERTIFIED TO ISO 13485. PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED.

Copyright © 2017 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, Scientists Helping Scientists, and MyeloCult are trademarks of STEMCELL Technologies Canada Inc. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.