

# Dissociation Reagents

## Collagenase B, ACF

Animal component-free collagenase for the digestion of native collagen fibrils

Catalog # 07438  
07439

100 mg  
1 g



Scientists Helping Scientists™ | [WWW.STEMCELL.COM](http://WWW.STEMCELL.COM)

TOLL FREE PHONE 1 800 667 0322 • PHONE +1 604 877 0713

[INFO@STEMCELL.COM](mailto:INFO@STEMCELL.COM) • [TECHSUPPORT@STEMCELL.COM](mailto:TECHSUPPORT@STEMCELL.COM)

FOR GLOBAL CONTACT DETAILS VISIT OUR WEBSITE

## Product Description

Collagenase B, Animal Component-Free (ACF) is obtained from cultures free of animal-derived materials. Collagenase is a protease consisting of a single polypeptide chain approximately 1,000 amino acid residues in length. Collagenase is capable of digesting native collagen fibrils commonly found in connective tissues and therefore is frequently used for tissue dissociation. Collagenase preparations contain the activity of several proteases, including collagenase, caseinase, clostripain, and trypsin (Kessler & Yaron). Collagenase B, ACF contains higher collagenase and caseinase levels than Collagenase A and is intended for use in applications where it is necessary to prevent the introduction of potential animal-derived pathogens.

## Product Information

Alternative Names:	Clostridium histolyticum collagenase; Collagenase B
Format:	Lyophilized powder
Storage:	Store at 2 - 8°C.
Stability:	Stable until expiry date (EXP) on label.
Reconstitution:	Dissociation reagents can be reconstituted in a balanced salt solution or buffer of choice.
Molecular Weight:	68 - 130 kDa
CAS Number:	9001-12-1
Optimum pH:	6.3 - 8.5
Cleavage Site:	-Pro-X-†-Gly-Pro-Y- : X = neutral Y = nonspecific

## Specifications

Source:	Clostridium histolyticum
Activity:	Collagenase: ≥ 300 CDU/mg dry weight (mgdw); Caseinase: ≥ 300 u/mgdw; Clostripain: ≤ 5.0 u/mgdw; Trypsin: ≤ 0.5 u/mgdw. See Notes for further information.

## Related Products

For a complete list of dissociation reagents, as well as related products available from STEMCELL Technologies, please visit our website at [www.stemcell.com](http://www.stemcell.com) or contact us at [techsupport@stemcell.com](mailto:techsupport@stemcell.com).

## Notes

### ACTIVITY UNITS

Collagenase: 1 collagenase digestion unit (CDU) equals 1  $\mu\text{mol}$  of L-leucine equivalents released from collagen in 5 hours at 37°C, pH 7.5.

Caseinase: 1 unit equals 1  $\mu\text{mol}$  of L-leucine equivalents released from 25 mg vitamin-free casein in 5 hours at 37°C, pH 7.5. Measures non-specific proteolytic activity.

Clostripain: 1 unit hydrolyzes 1  $\mu\text{mol}$  of N $\alpha$ -benzoyl-L-arginine ethyl ester (BAEE)/minute at 25°C at pH 7.6, after activation in 2.5 mM dithiothreitol (DTT).

Trypsin: 1 unit hydrolyzes 1  $\mu\text{mol}$  of BAEE/minute at 25°C at pH 7.6.

## References

- Bond MD & Van Wart HE. (1984) Characterization of the individual collagenases from *Clostridium histolyticum*. *Biochemistry* 23(13): 3085–91.
- Cherkashin AP et al. (2016) Expression of calcium-activated chloride channels Ano1 and Ano2 in mouse taste cells. *Pflugers Arch Eur J Physiol* 468(2): 305–19.
- Ghazanfari S et al. (2016) Modulation of collagen fiber orientation by strain-controlled enzymatic degradation. *Acta Biomater* 35: 118–26.
- Kempf H et al. (2015) Cardiac differentiation of human pluripotent stem cells in scalable suspension culture. *Nat Protoc* 10(9): 1345–61.
- Kessler E & Yaron A. (1973) A novel aminopeptidase from *clostridium histolyticum*. *Biochem Biophys Res Commun* 50(2): 405–12.
- Qi Z et al. (2016) TRPC3 regulates the automaticity of embryonic stem cell-derived cardiomyocytes. *Int J Cardiol* 203: 169–181.
- Sharifpanah F et al. (2015) Stimulation of vasculogenesis and leukopoiesis of embryonic stem cells by extracellular transfer RNA and ribosomal RNA. *Free Radic Biol Med* 89: 1203–17.
- Ugele B & Lange F. (2001) Isolation of endothelial cells from human placental microvessels: effect of different proteolytic enzymes on releasing endothelial cells from villous tissue. *In Vitro Cell Dev Biol Anim* 37(7): 408–13.
- Zhao T et al. (2016) Vascular endothelial growth factor-D mediates fibrogenic response in myofibroblasts. *Mol Cell Biochem* 413(1-2): 127–135.

STEMCELL TECHNOLOGIES INC.'S QUALITY MANAGEMENT SYSTEM IS CERTIFIED TO ISO 13485. PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED.

Copyright © 2016 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design and Scientists Helping Scientists are trademarks of STEMCELL Technologies Inc. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.