

ImmunoCult™ Human Treg Differentiation Supplement

Serum-free culture supplement for the differentiation of human naïve CD4+ T cells into regulatory T cells (Tregs)

Catalog # 10977 1 mL



Scientists Helping Scientists™ | WWW.STEMCELL.COM

TOLL FREE PHONE 1 800 667 0322 • PHONE +1 604 877 0713

INFO@STEMCELL.COM • TECHSUPPORT@STEMCELL.COM

FOR GLOBAL CONTACT DETAILS VISIT OUR WEBSITE

Product Description

ImmunoCult™ Human Treg Differentiation Supplement contains a human cytokine and small molecule formulated to promote the robust activation, expansion, and differentiation of peripheral blood-derived, naïve, CD4+ human T cells into regulatory T cells (Tregs).

ImmunoCult™ Human Treg Differentiation Supplement is intended for use with ImmunoCult™-XF T Cell Expansion Medium (Catalog #10981) and ImmunoCult™ Human CD3/CD28 T Cell Activator (Catalog #10971).

- Formulated for the induction of human Treg differentiation
- Optimized for use with ImmunoCult™-XF T Cell Expansion Medium
- Supports higher levels of memory T cells in comparison to competitor Treg differentiation kit
- Supports up to 20% higher proportions of CD4+CD25+FOXP3+T cells in comparison to competitor Treg differentiation kit
- Supplied as a 50X concentrate; after thawing and mixing, the tube contents can be added directly to medium

Properties

- Storage:** Store at -20°C.
- Shelf Life:** Stable for 12 months from date of manufacture (MFG) on label.
- Contains:**
- Recombinant human TGF-β1
 - All-trans retinoic acid

Handling / Directions For Use

PREPARATION OF COMPLETE Treg DIFFERENTIATION MEDIUM

Use sterile techniques to prepare complete Treg Differentiation Medium (ImmunoCult™ Human Treg Differentiation Supplement + ImmunoCult™-XF T Cell Expansion Medium). The following example is for preparing 50 mL of complete medium. If preparing other volumes, adjust accordingly.

1. Thaw ImmunoCult™ Human Treg Differentiation Supplement at room temperature (15 - 25°C) until just thawed. Mix thoroughly.
NOTE: If necessary, centrifuge for 30 seconds to recover liquid from cap.
NOTE: If not used immediately, store supplement at 2 - 8°C for up to 1 month. Do not re-freeze.
2. Add 1 mL of ImmunoCult™ Human Treg Differentiation Supplement to 49 mL of ImmunoCult™-XF T Cell Expansion Medium (Catalog #10981). Mix thoroughly.

DIFFERENTIATION TO Tregs

1. Start with purified human naïve CD4+ T cells at 1×10^6 cells/mL in complete Treg Differentiation Medium.
2. Activate human naïve CD4+ T cells using ImmunoCult™ Human CD3/CD28 T Cell Activator (Catalog #10971) or other desired reagents to activate T cells. Incubate cells at 37°C and 5% CO₂.
3. Adjust the cell density to 1×10^6 cells/mL every 3 - 4 days as needed with the addition of fresh complete Treg Differentiation Medium for 7 days.
4. On Day 7, differentiated Tregs are ready to be used in the desired application.

STEMCELL TECHNOLOGIES INC.'S QUALITY MANAGEMENT SYSTEM IS CERTIFIED TO ISO 13485. PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED.

Copyright © 2017 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, Scientists Helping Scientists, and ImmunoCult are trademarks of STEMCELL Technologies Canada Inc. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.