

STEMdiff™ Cardiomyocyte Support Medium

Medium for thawing and culturing hPSC-derived cardiomyocytes



Scientists Helping Scientists™ | www.stemcell.com

TOLL FREE PHONE 1 800 667 0322 • PHONE +1 604 877 0713

INFO@STEMCELL.COM • TECHSUPPORT@STEMCELL.COM

FOR GLOBAL CONTACT DETAILS VISIT OUR WEBSITE

Catalog #05027

250 mL

Product Description

STEMdiff™ Cardiomyocyte Support Medium is a versatile medium that limits stress on human pluripotent stem cell (hPSC)-derived cardiomyocytes when transitioning from cryopreservation to thawing and from harvesting to replating of hPSC-derived cardiomyocytes. After thawing or replating, the functional capacity of hPSC-derived cardiomyocytes is retained and cells can be used in various downstream applications and analyses. The hPSC-derived cardiomyocytes can be further maintained long-term using STEMdiff™ Cardiomyocyte Maintenance Kit (Catalog #05020).

Product Information

PRODUCT NAME	CATALOG #	SIZE	STORAGE	SHELF LIFE
STEMdiff™ Cardiomyocyte Support Medium	05027	250 mL	Store at -20°C.	Stable for 12 months from date of manufacture (MFG) on label.

Materials Required But Not Included

PRODUCT NAME	CATALOG #
Corning® Matrigel® hESC-Qualified Matrix	Corning 354277
STEMdiff™ Cardiomyocyte Maintenance Kit <ul style="list-style-type: none">STEMdiff™ Cardiomyocyte Maintenance Basal MediumSTEMdiff™ Cardiomyocyte Maintenance Supplement (50X)	05020
Trypan Blue	07050

Directions for Use

Please read the entire protocol before proceeding. Use sterile techniques when performing the protocols below.

A. THAWING hPSC-DERIVED CARDIOMYOCYTES

Frozen hPSC-derived cardiomyocytes should be thawed and plated onto Corning® Matrigel®-coated cultureware. For coating plates with Corning® Matrigel®, refer to the Technical Manual: Maintenance of Human Pluripotent Stem Cells in mTeSR™1 (Document #28315) or TeSR™-E8™ (Document #DX20809) available at www.stemcell.com or contact us to request a copy.

For storage, stability, and preparation instructions for STEMdiff™ Cardiomyocyte Maintenance Medium, refer to the corresponding Product Information Sheet (Document #DX21496), available at www.stemcell.com or contact us to request a copy.

- Coat a 12-well tissue culture plate with Corning® Matrigel® hESC-Qualified Matrix and bring to room temperature (15 - 25°C) for at least 1 hour prior to use.
- Thaw STEMdiff™ Cardiomyocyte Support Medium at room temperature (15 - 25°C) or overnight at 2 - 8°C. Warm to room temperature.
- Thaw hPSC-derived cardiomyocytes in a 37°C water bath by gently shaking the cryovial continuously until only a small frozen cell pellet remains.
- Add 5 - 7 mL of STEMdiff™ Cardiomyocyte Support Medium to a 15 mL conical tube (e.g. Catalog #38009).
- Using a 2 mL pipette, gently transfer the contents of the cryovial to the tube from step 4.
- Centrifuge the cells at 300 x g for 5 minutes at room temperature (15 - 25°C).

7. Aspirate the supernatant and gently add 1 - 2 mL of STEMdiff™ Cardiomyocyte Support Medium to resuspend cells.
8. Perform a cell count using Trypan Blue and a hemocytometer.

B. CULTURING hPSC-DERIVED CARDIOMYOCYTES

1. Aspirate Corning® Matrigel® from a pre-coated 12-well tissue culture plate (from section A step 1). Add 2 mL of STEMdiff™ Cardiomyocyte Support Medium per well.
2. Add cells at a density appropriate for downstream assays or other applications. Incubate at 37°C for 24 hours.
3. Remove medium and add 2 mL of STEMdiff™ Cardiomyocyte Maintenance Medium per well. Incubate at 37°C.
4. Every 2 days, perform a full medium change with 2 mL of STEMdiff™ Cardiomyocyte Maintenance Medium per well.

STEMCELL TECHNOLOGIES INC.'S QUALITY MANAGEMENT SYSTEM IS CERTIFIED TO ISO 13485. PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED.

Copyright © 2019 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, Scientists Helping Scientists, and STEMdiff are trademarks of STEMCELL Technologies Canada Inc. mTeSR, TeSR, and E8 are trademarks of WARF. Corning and Matrigel are registered trademarks of Corning Incorporated. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.